



668 N. Coast Hwy. Suite 1177 Laguna Beach, CA. 92651  
P. 1-888-402-2778 F. 1-888-901-4166  
[www.m2associates.com](http://www.m2associates.com) [tom@m2associates.com](mailto:tom@m2associates.com)

## Olympus BH2 RLA Vertical Illuminator Operation



Push Pull Rod – Push in for Dark Field Observation. Pull out for Bright Field Operation.  
Note: Some RLA's have the ND Filter removed. If it is present, simply press down on the ND Filter rod and it will move out of the way. Olympus only put this on the RLA illuminator.

Aperture Diaphragm – Adjust this to the Numerical Number of the Objective Lens. For example, if you are using a 100X objective the Numerical Aperture is .90. To optimize the resolution of the lens close, the Aperture down 90%. You can visually see resolution improve as you close this Aperture down.